

MAGPIX® Drive Fluid 4 Pack

Product Information Sheet



Summary

Use MAGPIX® Drive Fluid as the delivery and flush medium, which carries the sample to the optics component of the MAGPIX based instruments.

Safety Precautions

Avoid contact with skin and eyes. A Safety Data Sheet (SDS) is available upon request.

Ingredients

MAGPIX® Drive Fluid contains sodium chloride, sodium phosphate, and an antimicrobial in water.

Product Storage

Store at 15°C to 30°C. Do not freeze.

Limitations

You must follow these product information sheet instructions. Reliability of results cannot be guaranteed if you deviate from these instructions. When stored at 15°C to 30°C, the product should perform as expected up to the expiration date stated on the container label.

Procedure

1. Open the carton of drive fluid. Open the MAGPIX® instrument fluid compartment and remove the empty drive fluid bottle and dispose of properly.
2. Slide the new drive fluid bottle part in, then firmly grip the seal tab while holding the bottle and remove the foil seal.
3. Extend the drive fluid tubing from the fluid compartment and carefully insert the plug in the drive fluid bottle opening ensuring the tubing is pointing straight down into the bottle indentation.
4. Slide the drive fluid bottle in place inside the compartment. The bottle should fit into the indents of the tray. Check that the waste container is empty.
5. Prime the unit per the user manual's instructions.

NOTE: Product Information Sheets are available in other languages upon request. Contact Luminex Technical Support for information or access on website at <http://www.luminexcorp.com>. Then, using the Search feature, search for the desired product information sheet.

Refer to the user manual for proper use of the MAGPIX Drive Fluid as applicable.

RUO For Research Use Only. Not for use in diagnostic procedures.
89-60000-00-075 Rev C
January 2020

REF Product No. MPXDF-4PK
Quantity: 4 x 700 mL

To order more product:
Email: orders@luminexcorp.com
Fax: 512-219-0544

© 2010-2020 Luminex Corporation

Technical Support

Telephone: 512-381-4397
North America Toll Free: 1-877-785-2323
International Toll Free: + 800-2939-4959
Email: support@luminexcorp.com
www.luminexcorp.com



Luminex Corporation

12212 Technology Blvd.
Austin, TX 78727
U.S.A.